## S90322 REDI-SEP®

For use in emergency  
eye wash stations

A preservative  
for potable water

8 Fluid Ounces (236 ml)

### Product information

Bottle Contents: Chlorhexidine Gluconate 1% w/v, CAS 18472-51-0, Propylene Glycol CAS 57-55-6

Uses: An additive/conditioner combined with potable water, which extends the water change cycle on portable eye wash stations.

Warnings: Do not use if seal is broken or missing. Do not allow contact with eyes prior to mixing with potable water according to the directions. In case of contact with eyes in undiluted form, flush with clean water. If you experience any eye pain, change in vision, continued redness or irritation, or if condition worsens after treatment at the eye wash station, immediately consult a physician.

Keep this and all chemicals out of reach of children. Do not change the dilution or add other ingredients.

Directions: Clean portable, self-contained station with potable water and rinse. Partially fill station with potable water. Add entire contents of 8 oz. REDI-SEP® bottle. Finish filling tank according to the filling instructions supplied with eye wash. Date and initial inspection tag.

REDI-SEP® when mixed as directed, preserves eye wash station water volumes at minimum levels of 4 gallons (15 liters) and max levels of 20 gallons (75 liters). Eye wash station lines should be flushed weekly with a 15 second flush to ensure proper operation.

All test water should be replaced with clean potable water. Replacement potable water for these tests will not reduce the effectiveness of the REDI-SEP® to the station. Refer to instructions supplied with eye wash for further testing procedures.

It is recommended that the station be cleaned and the water replaced at a minimum of every six months. Examination at weekly intervals which give evidence of any particulates or discoloration suggests the water be replaced immediately.

Other Information: Store at room temperature 15°-30° C (60°-80° F)

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